

# Rely+On™ Virkon™


Broad Spectrum  
Virucidal Disinfectant

## How to prepare and use a Rely+On™ Virkon™ disinfectant solution

### Preparing a 500ml 1:100 disinfectant solution with Rely+On™ Virkon™ 5g tablets


#### STEP 1

Wear appropriate personal protective equipment before preparation and use of disinfectant solutions

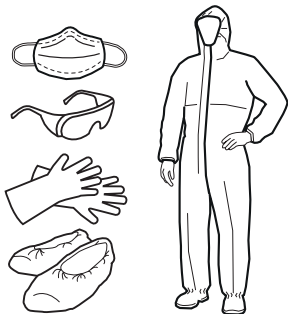

#### STEP 2

Fill a 500ml trigger spray bottle with warm water

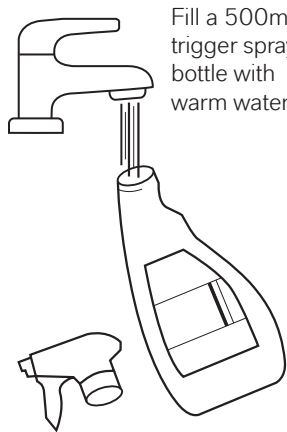


#### STEP 3

Add 1 x 5g Rely+On™ Virkon™ tablet and wait for it to dissolve completely



#### STEP 4


The Rely+On™ Virkon™ disinfectant is now ready for use

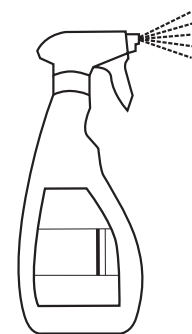

### Preparing a 500ml 1:100 disinfectant solution with Rely+On™ Virkon™ powder concentrate


#### STEP 1

Wear appropriate personal protective equipment before preparation and use of disinfectant solutions

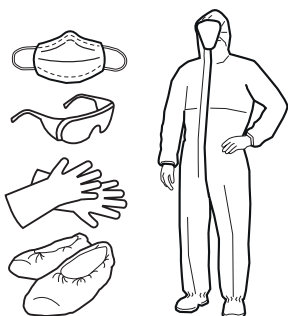


#### STEP 2

Fill a 500ml trigger spray bottle with warm water

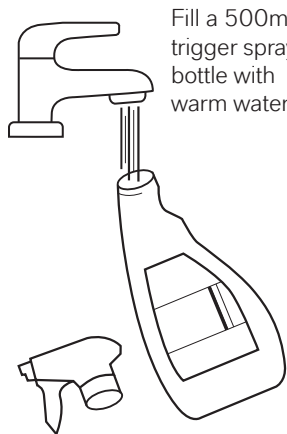

#### STEP 3


Weigh out 5g Rely+On™ Virkon™ powder concentrate

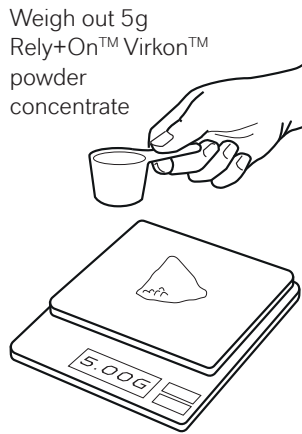

#### STEP 4

Add the 5g **+5g** to the bottle and wait for it to dissolve completely


The Rely+On™ Virkon™ disinfectant is now ready for use

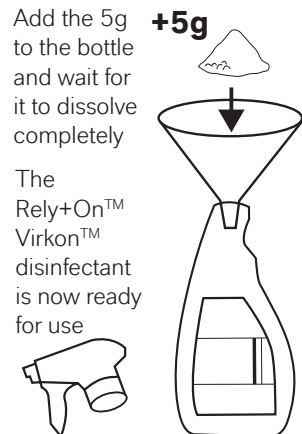

### ALTERNATIVE METHOD USING MEASURING JUG:


#### STEP 1

Fill a jug with 500ml warm water

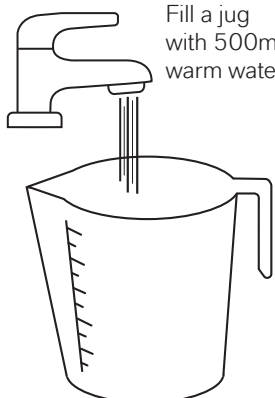


#### STEP 2

Add 5g Rely+On™ Virkon™ powder concentrate to jug, and stir to dissolve completely

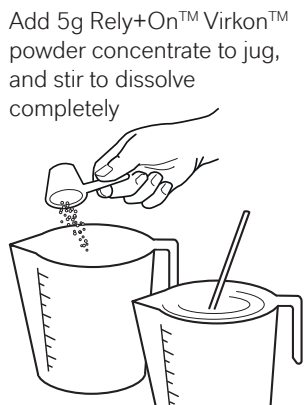


#### STEP 3

Leave for 5 minutes, then pour solution into spray bottle

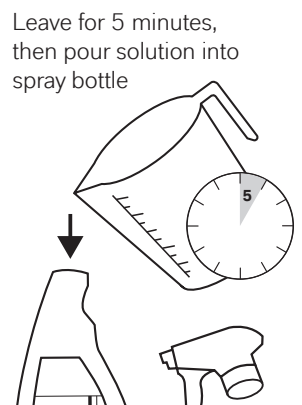

#### STEP 4

The Rely+On™ Virkon™ disinfectant is now ready for use

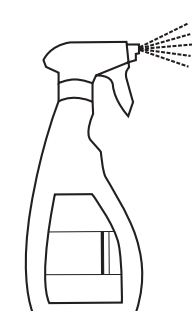

## Applying Rely+On™ Virkon™ disinfectant solution via a trigger spray bottle


### STEP 1

Spray Rely+On™ Virkon™ solution onto the surface to be disinfected



### STEP 2

Wait a minimum of 10 minutes†

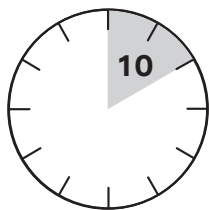


### STEP 3

Spray all surfaces that have had disinfectant applied with fresh clean water



### STEP 4

Then wipe surface dry with a clean cloth

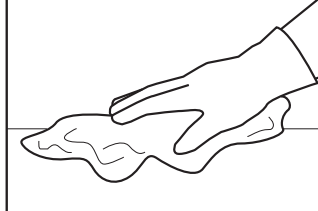


† 1 minute for control of SARS-COV-2 virus

## Preparing 1:100 disinfectant solutions with Rely+On™ Virkon™ powder concentrate


### STEP 1

Wear appropriate personal protective equipment before preparation and use of disinfectant solutions

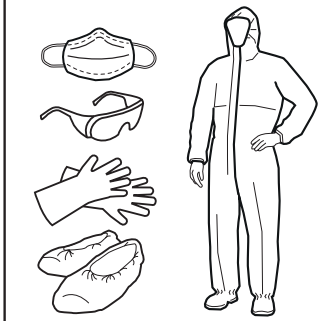


### STEP 2

Fill a suitable container with the required volume (e.g. 5 litres) of warm water

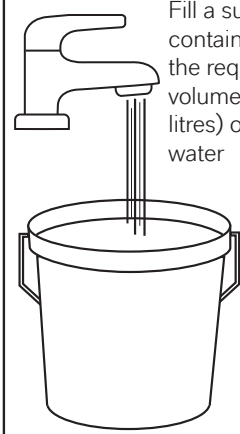


### STEP 3

Weigh out 10g of Rely+On™ Virkon™ powder concentrate for every 1 litre of water

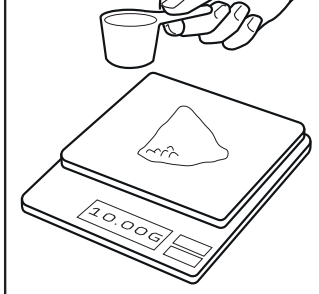

### STEP 4


Add the Rely+On™ Virkon™ powder to the water and then stir until completely dissolved. The disinfectant solution is now ready to use

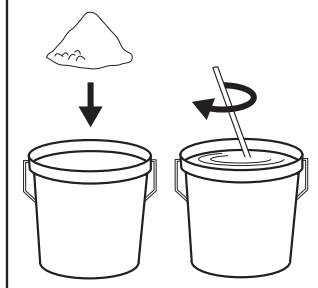

## Applying Rely+On™ Virkon™ disinfectant solutions


### STEP 1

To disinfect large areas of floor space, apply Rely+On™ Virkon™ solution using a mop and bucket

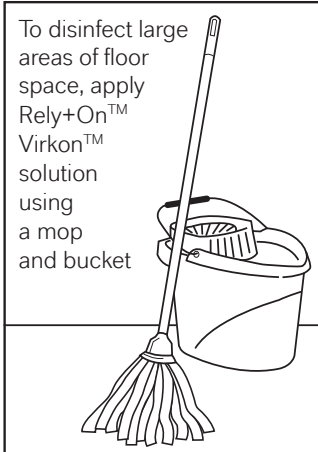


### STEP 2

To disinfect walls, doors, and other large surfaces, use a bucket and sponge or cloth to apply Rely+On™ Virkon™ disinfectant solution

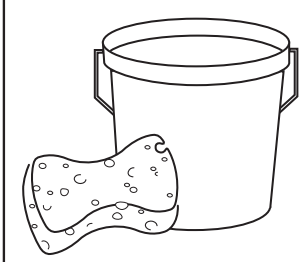

### STEP 3

Wait for a minimum of 10 minutes†

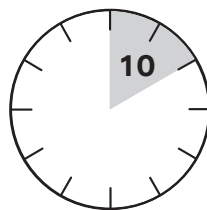

### STEP 4

Rinse all surfaces with clean water when materials compatibility is of concern



† 1 minute for control of SARS-COV-2 virus

\* Not for use for the disinfection of medical devices.

\*\* Users should always check for compatibility with any sensitive materials before use. Do not use on any soft metals such as brass or copper. Not recommended for use on soft furnishings, textiles or carpets.

The specified uses and registered claims may vary from country to country. Please contact LANXESS to verify country-specific approved uses.

**Use biocides safely. Always read the label and product information before use.**



Antec International Limited  
Windham Road, Chilton Industrial Estate  
Sudbury, Suffolk CO10 2XD,  
United Kingdom

Tel: +44(0)1787 377305  
relyondisinfection@lanxess.com  
www.relyondisinfection.com

©2020 LANXESS. Rely+On™ Virkon™, LANXESS™ and the LANXESS Logo are trademarks of LANXESS Deutschland GmbH or its affiliates. All trademarks are registered in many countries worldwide.

How to prepare/Rely+On™Virkon™/UK/01.10.20/V6/G